

SCAQMD M313-91

VOC by GC/MS/FID

Technical Discussion
September 5th, 2013

1

Agenda

Method Endpoint
& Other Policy Decisions

Following
Meetings

2

Agenda

Introductions

Technical Presentation 30 minutes

Comparative Data 20 minutes

Analytical Concerns
(Previous Meeting) 15 minutes

Take Comments 30 minutes

Extend Meeting? Set Next Agenda?

3

Method Context

M24
Non-Volatiles
Density
Karl-Fischer
Exempts (if necessary)

Method 24 Analysis Completed

Every Sample

4

M313 Usage

< 150 g/L Sample Concentration

Inconsistent M24 Weighings

District Rules Or Protocols (CAS)

Limited Applicability

5

Method Procedure

Calibration

Sample Pre-Screening

Sample & QC Preparation

Sample Processing

6

Calibration

Calibrated Compound	RRF
Ethylene Glycol	0.5
Propylene Glycol	0.7
Texanol	1.4
1-Butanol	1.3

Pre-Calibrate
Commonly Seen
Compounds

Iterative In Nature

7

Sample Pre-Screening

Required Due To Pre-Calibration Requirement

Scenario #1

Sample
Pre-screened



No New
Compounds
Detected

Proceed To Sample Analysis

8

Sample Pre-Screening

Required Due To Pre-Calibration Requirement

Scenario #2

Sample
Pre-screened



At Least 1 New
Compound
Detected

Must Pre-Calibrate New Compounds

9

Calibration

Calibrated Compound	RRF
Ethylene Glycol	0.5
Propylene Glycol	0.7
Texanol	1.4
1-Butanol	1.3
Decane	2.0

Scenario #2

New Compounds
Calibrated & Added To
Calibration List

Proceed To Sample Analysis

10

Sample Analysis



Clear Lacquer

~ 30 Grams



Lacquer Aliquot

11

Sample Analysis



Lacquer Aliquot



IPA
Heptane
Triglyme
DIIBA



Spiked Lacquer
Mixed Well

12

Surrogate Spikes

Compound	Characteristic	Retention	Functional Group
IPA	Small, Polar	Early	Alcohol
Heptane	Small, Non-Polar	Early-Mid	Alkane
Triglyme	Large, Polar	Mid-Late	Poly-ether
DIIBA	Large, Non-Polar	Late	Ester

13

Sample Analysis



Spiked Lacquer
Mixed Well




2.5 Grams
Of Sample
& EGDE
(Internal Standard)



Diluted Lacquer
To 25 mL
With Solvent

14

Surrogate Addition

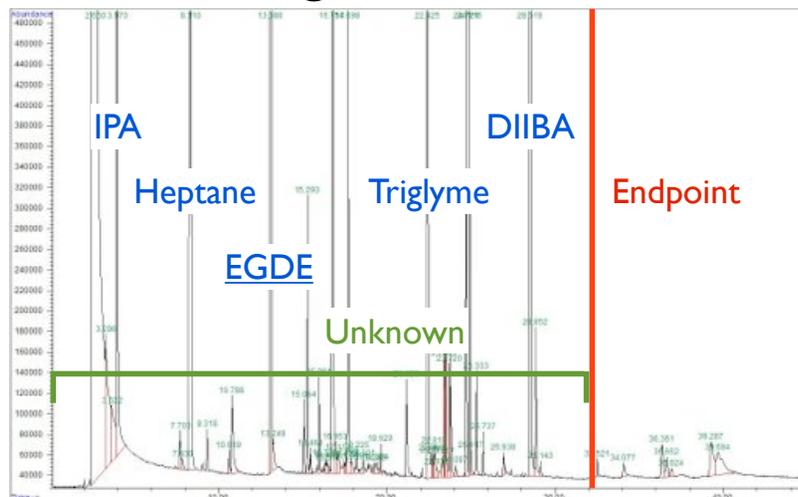

15

Batch Standards



Run With Every Sample Set

Reagent Blank (RB)

Continuing Spike Verification (CSV)

Target Compound Mix (TCM)


Instrument Optimization Mix (IOM)

16

Instrument Optimization Mix

Component Evaluates

C6 - C15	Discrimination
BFB	Fragmentation
Methyl Palmitate	Endpoint
Ethylene Glycol	Resolution
Propylene Glycol	Resolution
Triglyme (0.1 g/L)	LOD

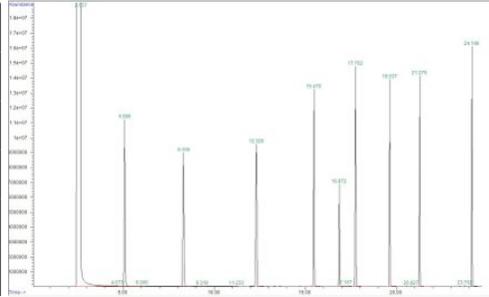

17

Processing

Contaminant Subtraction


Relative Response Factor Assignment

Sum VOC Totals

18

Triglyme Default Compound

Every Peak First Quantified As

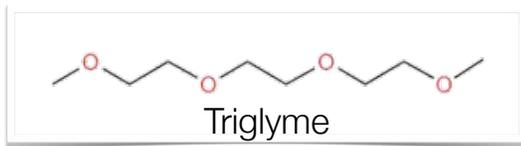

Concentration "As Triglyme"
Dictates Next Step

19

Surrogate Compound Usage

Concentration
As Triglyme

< 0.1 g/L

Action

Ignore Peak

20

Surrogate Compound Usage

**Concentration
As Triglyme**

0.1 - 1 g/L

Action

Use "As Triglyme" Concentration
For Final VOC Tally

21

Surrogate Compound Usage

**Concentration
As Triglyme**

1 - 2 g/L

Action

Calculate Using RRF Of Identified Compound
Or A Similar Compound
[Same Functional Groups; 1 Carbon Difference]

22

Surrogate Compound Usage

**Concentration
As Triglyme**

2+ g/L

Action

Calculate Using RRF Of
Identified Compound Only

23

QC Recoveries

24

Surrogate Spikes

Average % Recovery

	D 6886	Intralab
EGDE (IS)	94	92
IPA	101	100
Heptane	94	99
Triglyme	98	99
DIIBA	99	93

25

M24 Comparative Data

ASTM D6886 Interlab
2010

26

	M24 (g/L)	M313 (g/L)
#1	-22	6
#2	14	3
#3	-15	13
#4	-34	5
#5	-12	17
#6	-25	5

“Near 0 g/L”
As M24 (Coating)

ASTM D6886
Interlab 2010

27

	M24 (g/L)	M313 (g/L)
#1	35	70
#2	-6	51
#3	-7	53
#4	-8	55
#5	-7	55
#6		

“50 g/L”
As M24 (Coating)

ASTM D6886
Interlab 2010

28

M24 Comparative Data

Random Samples
2008 - 2012

(g/L)

	M24	M313	Type
#1	5	6	Latex Ceiling Paint
#2	13	39	Acrylic Sealer
#3	16	13	Enamel Undercoater
#4	25	13	Roof Coating
#5	25	21	Traffic Paint
#6	33	44	Acrylic Semi-Gloss
#7	36	36	Matte Finish
#8	42	41	Concrete Coating
#9	47	83	Polyurethane Sealer
#10	53	103	Flat Varnish

Sample

0 - 50 g/L
Samples As
M24 (Material)

Random Samples
2008 - 2012

(g/L)

	M24	M313	Type
#1	55	41	Concrete Coating
#2	59	46	Exterior Latex
#3	60	77	High Gloss Enamel
#4	63	43	Wood Watersealer
#5	72	57	Traffic Paint
#6	75	74	Penetrating Sealer
#7	78	48	Sealer & Finish
#8	81	26	Traffic Paint
#9	82	81	Flat Varnish

Sample

50 - 100 g/L
Samples As
M24 (Material)

Random Samples
2008 - 2012


(g/L)

	M24	M313	Type
#1	106	127	Concrete Sealer
#2	110	90	Exterior House Stain
#3	125	80	Flat Varnish
#4	136	110	Traffic Paint
#5	143	123	Clear Wood Finish

Sample

100 - 150 g/L
Samples As
M24 (Material)


Random Samples
2008 - 2012

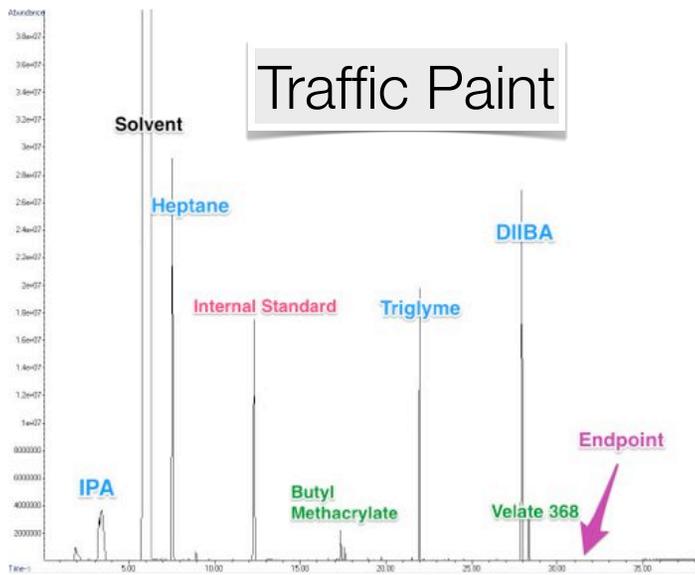

Traffic Paint

M24
81 g/L

M313
26 g/L

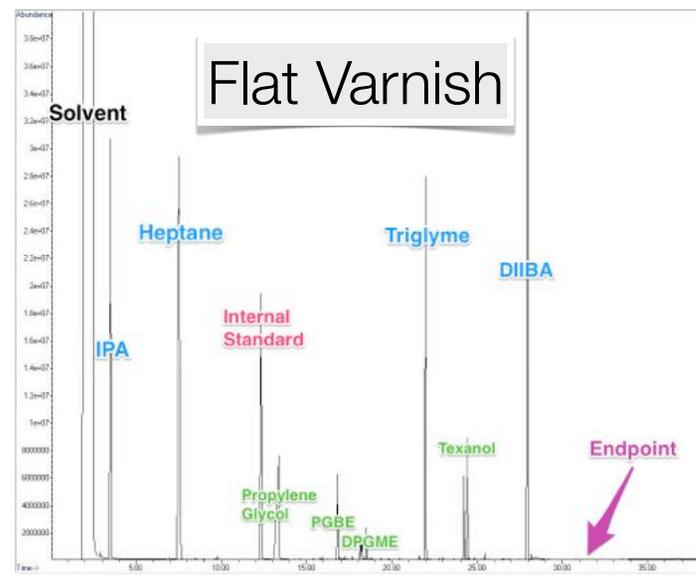

Flat Varnish

M24
53 g/L

M313
103 g/L

Analytical Concerns

Inlet Breakdown

Inlet Cooling With Reference Compound

New Liner Every Sequence

DIIBA Reactivity

Formulation Data

GC/MS Alternative

FID-Only Alternative Method

“Known” Formulations

Reduced QC; IOM mandatory

Will Discuss With EPA

37

What about D 6886?

Potentially Acceptable Substitute

Endpoint Differences

Use Triglyme Default Compound; IOM Mandatory

Will Discuss With EPA

38



M313 Test Methodology

Joan Niertit, Principal Chemist
(909) 396-2174
jniertit@aqmd.gov

Brad Parrack, Air Quality Chemist
(909) 396-3071
bparrack@aqmd.gov

39